

PHYWE Systeme GmbH & Co. KG
Robert-Bosch-Breite 10
37079 Göttingen
Germany

Telefon +49 (0) 551 604-0
Fax +49 (0) 551 604-107
E-mail info@phywe.de
Internet www.phywe.com

Operating instructions



Fig. 1: PHYWE Demo multimeter ADM 3 light, 13830-00

TABLE OF CONTENTS

- 1 SAFETY PRECAUTIONS
- 2 PURPOSE AND INTENDED USE
- 3 FUNCTIONAL AND OPERATING ELEMENTS
- 4 OPERATING INSTRUCTIONS
- 5 HANDLING
- 6 FIRMWARE UPDATE
- 7 TECHNICAL DATA
- 8 SCOPE OF SUPPLY
- 9 ACCESSORIES
- 10 REPLACING THE OVERLOAD SAFETY DEVICE
- 11 WARRANTY
- 12 DISPOSAL
- 13 CLEANING
- 14 EXPLANATION OF THE SYMBOLS

1 SAFETY PRECAUTIONS



on!Attention

The instrument complies with the safety requirements of DIN EN 61010-1. The instrument can be used safely and without risk if the following safety instructions are strictly adhered to.

- Read the operating instructions thoroughly and completely prior to starting the instrument. This is important for your own protection and for avoiding damage to the instrument.
- In the event of a change of ownership, these operating instructions must be handed over with the instrument.
- Use the instrument only for its intended purpose. Do not exceed the maximum values stated in section 7 „Technical data“.
- If the instrument is switched off by manual selection of the "OFF" mode or if it is switched off automatically due to standby or for deep-discharge protection, the measuring leads must be disconnected.
- The instrument is solely intended for use in dry rooms where there is no risk of explosion.
- Position the instrument on a table in an upright position for using or charging it.
- The hollow grip is solely intended for transporting the switched-off instrument to its location of use. Do not swing the instrument around when transporting it.
- Do not open the instrument. Only PHYWE is authorised to change the batteries.
- Only use the approved accessories.
- Do not connect any devices to the instrument other than the ones that are intended for this purpose.

- Prior to opening the back panel, it is mandatory to switch the instrument off and to disconnect all of the measuring leads.
- Do not use the instrument unless its housing is closed and undamaged.
- When measuring voltages above 30 V, only use connecting leads with safety connectors (safety connecting leads with fixed insulating sleeves).
- Ensure that no liquids or objects penetrate the instrument through its slots.
- Protect the instrument against shocks, dust, moisture vapours, direct exposure to sunlight and extremely high temperatures.
- Clean the instrument only with a slightly moist, lint-free cloth. Do not use any aggressive cleaning agents or solvents.
- Do not use the instrument if you notice signs of damage to the instrument, power supply unit, mains power cable or measuring leads. Do not use the instrument if it does not work correctly.

2 PURPOSE AND INTENDED USE

The PHYWE Demo multimeter ADM 3 light is a modern, universal multi-range meter that has been specifically developed for teaching purposes. The integrated measuring amplifier enables voltage and current measurements over an extremely wide range.

The instrument is characterised by a high input resistance in terms of the voltage measuring ranges and a particularly small voltage drop in terms of the current measuring ranges. It has eight selectable scales, one of them with central zero. In addition, all of the measurements can also be displayed as digital numbers.

The instrument is designed for measuring applications in test and measuring circuits that are connected to the mains power supply directly or indirectly. For test and measuring circuits, that are connected to the mains power supply directly, the instrument is designed for measurement category II 300 V (abbrev. CAT II 300 V). Allowed are measurements in circuits that are connected via a single-phase power plug to the low voltage system installation (incl. overcurrent protective element).



Attention!

Do not perform measurements of the building wiring installation. (e.g. fuse box, electric meter, service box, low voltage overhead lines)!

3 FUNCTIONAL AND OPERATING ELEMENTS

Fig. 2 shows the functional and operating elements at the front of the instrument and Fig. 3 those at the back.

3.1 Front

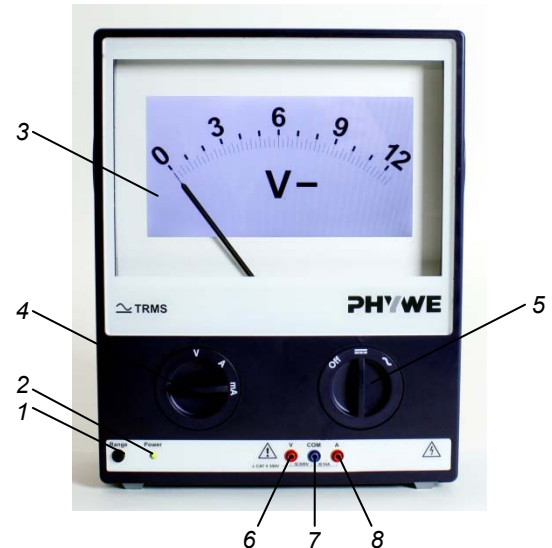


Fig. 2: Functional and operating elements, front

- 1 **"Range" scale button**
for selecting scales and digital display.
- 2 **"Power" LED**
for indicating the operating state:
On: Ready
Flashing: Standby
Off: Rotary switch set to OFF or protection against deep discharge
- 3 **Display**
for displaying the measured values in an analogue or digital form.
- 4 **Measuring range selector**
for selecting the measuring range:
Voltage V – current A / mA
- 5 **Operating mode selector**
for switching the device off (OFF) and for selecting the type of current (--- (direct current DC) or ~ (alternating current AC)).
In the case of AC, true root mean square (TRMS) values are measured. The input signal thereby is not limited to sinusoidal waves but measures the correct root mean square value for any given waveform. Also alternating signals with a share of direct signals are included (AC+DC measurement).
- 6 **4 mm safety socket V (red)**
for V measurements.
In case of DC voltage, connect the positive terminal here.
- 7 **4 mm safety socket COM (blue)**
reference point for the measuring sockets V and A.
- 8 **4 mm safety socket A (red)**
for all current measuring ranges.
In case of DC current, connect the positive terminal here.

3.2 Back

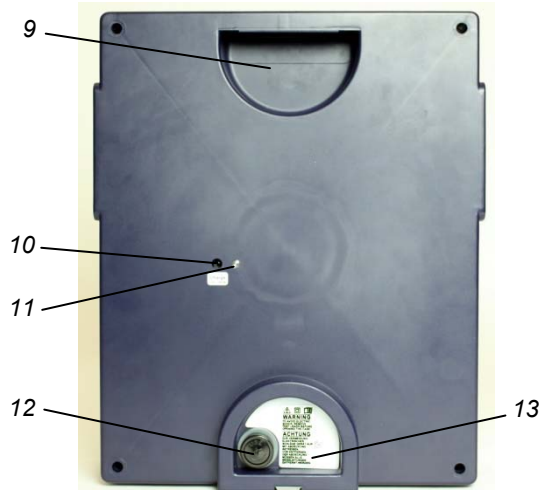


Fig. 3: Functional and operating elements, back

- 9 *Hollow grip*
for transportation.

- 10 *Charging connector*
for connecting the supplied charger.



Attention!

Only use the charger stated in the chapter 8 "Scope of supply" for charging the integrated batteries.

- 11 *Charge status LED*
for indicating the charge status of the battery:
yellow/orange: Charging
green: Fully charged (see sec. 4.2 "Charging")

- 12 *Fuse holder*
for overload protection.



Attention!

Do not use the instrument unless the fuse holder is closed!

- 13 *Cover*



Attention!

Do not use the instrument unless the cover is closed and fully intact!

Prior to removing the cover, switch the instrument off and disconnect all of the measuring leads. (see section 6.1 "Removing the cover")

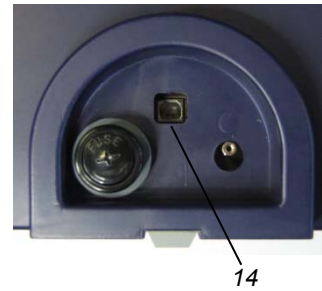


Fig. 4: Functional elements below the cover

14 *USB port*

for connecting a PC for a firmware update with the PHYWE ADM 3 Update Program. (see section 6.2 "Firmware update")

4 OPERATING INSTRUCTIONS

The instrument fulfils the technical requirements that are summarised in the current directives of the European Community.

The connected cables must not be longer than 2 m. Only the two sockets that are required for a particular measurement unit must be used.

4.1 Area of application

Near strong high-frequency transmitters, e.g. radio sets, the operating behaviour of the instrument may exceed the stated specifications.

The instrument must be used under the supervision of an expert and in an electromagnetically controlled environment at research, teaching and training facilities (schools, universities, institutes and laboratories).

This means that, in such an environment, radio transmission devices, e.g. mobile phones, should not be used in the direct vicinity of the instrument.

Electrostatic charges or similar electromagnetic phenomena (HF, bursts, indirect lightning discharge, etc.) may affect the instrument so that it will not work within the specified data range.

The following measures reduce or eliminate any potential interference:

Avoid carpets; provide equipotential bonding; perform the experiments on a conductive, earthed surface, use shields and shielded cables. Do not use radiofrequency transmitters (radio sets, mobile phones) in the direct vicinity of the instrument.

After a complete shut-down of the instrument, switch it off and then on again by way of the operating mode selector (5).

4.2 Charging

The instrument can be charged when it is switched on or off. During charging, the following symbol will be displayed:



To ensure trouble-free operation and a long service life of the instrument, the battery must be charged regularly. The operating time in the battery mode depends on the ambient temperature and selected measuring ranges.

During loading, interference signals are generated which superimpose the measurement signal. It is therefore recommended, especially with small measuring ranges, to measure exclusively in battery mode.

5 HANDLING

5.1 Start-up

To switch the instrument on, select the desired type of current with the aid of the operating mode selector (5). It will be ready for operation after a turn-on period of several seconds. The measuring range selector (4) and scale button (1) are used to select the measurement unit and scale and, thereby, the desired measuring range. After this step, the measuring circuit can be connected.

5.2 Buzzer

The buzzer is used for signalling the following operating states:

- **Short beep:** Overrange, >5-10 % beyond the scale range.

Display:



Elimination of the problem:

The overrange state must be eliminated by changing the measuring range (scale) or by reducing the measured variable!

- **Continuous sound:** overrange and deactivation of the measuring amplifier (to protect the device). The pointer returns to "0".

Display:



Elimination of the problem:

The overrange state must be eliminated by reducing the measured variable!

After a waiting period of several minutes, the measuring amplifier can be switched on again by changing the measuring range (selector 4) or operating mode (selector 5).

5.3 Standby

To increase the service life of the battery, the instrument is equipped with an automatic shut-off system that switches the instrument off approx. 60 minutes after the last intervention by an operator. The fact that the instrument has been switched off is indicated by the inactive display and a flashing "Power" LED (2).

In order to reactivate the instrument, the scale button (1), the measuring range selector (4) or the operating mode selector (5) must be actuated.

To switch the instrument completely off, the operating mode selector must be set to "OFF".

5.4 Battery mode

The instrument can be operated with a connected charger or by way of the integrated battery.

When there is only approx. 1 hour of battery life left, the following symbol will be displayed:



When the batteries are discharged, the instrument switches itself off so that there will be no deep discharge and, thereby, no damage to the batteries. After the automatic shut-off, the instrument should be switched off with the aid of the operating mode selector (5) (OFF).

The instrument can be charged for as long as desired. Once the battery is fully charged, the charge will be maintained.

5.5 Correct measurement of unsmoothed, pulsating DC voltages and currents

When measuring pulsating DC voltages and currents like the ones obtained from normal variable transformers with rectifiers, moving coil instruments are subject to a fundamental error of measurement: These instruments indicate the mean value instead of the root-mean-square (rms) value. In the case of a pulsating DC voltage obtained by way of a bridge rectifier, this mean value is almost exactly 11 % lower than the rms value.

However, the instrument still enables the exact measurement of the rms value of pulsating DC voltages and currents. Simply switch the instrument to the AC mode . In the case of sinusoidal pulsating measurement signals, the instrument indicates the exact rms value. The exact rms value is also obtained in the case of distorted sinusoidal signals. (For pure DC currents, the operating mode selector (5) should always be set to since, otherwise, the measurement accuracy will be reduced.)

5.6 Measurement of non-sinusoidal AC voltages and currents

To measure AC voltages and currents, the voltages and currents are rectified in the ADM 3 light in the mode by way of a true RMS-to-DC converter. The advantage of this method is that, even in the case of signals that deviate substantially from the sinusoidal form, exact rms values will be displayed. For signals with a crest factor (this is the ratio of peak value to RMS value) of 4, hardly any influence of the waveform on the measurement result can be observed.

5.7 Overload protection

The instrument is equipped with numerous protective devices to protect it against damage caused nearly all conceivable types of faults.

Voltage overload

The "V" socket (6) is protected against mains voltage in all of the measuring ranges and protected by a surge protector against higher voltages (short-time up to 1kV). For non-hazardous voltages that are safe to touch, the protective function is effective up to 10 kV.

A voltage is considered to be non-hazardous (i.e. safe to touch) if it complies with the following conditions (according to standard DIN EN 61010-1):

- the values are smaller than 33 V_{rms} in the case of AC voltage or
- smaller than 70 V in the case of DC voltage or
- the current (terminated with 2 kΩ) is smaller than 2 mA in the case of direct current or
- the capacitive charge is smaller than 45 μC (valid up to U<15 kV) or
- the stored energy value is smaller than 350 mJ (valid for U>15 kV).

Current overload

The socket „A“ (8) is protected against current overload by two protective measures: An overload cut-off (thermal cut-off) and a fuse for high currents. Due to overrange, the measuring circuit "A" will be cut-off not later than 3 s and the measuring circuit "mA" not later than 15 s. Depending on the overload of the current measuring ranges, a thermo switch may temporarily disrupt the circuit.

After a cooling period of several minutes, the current measurement function can be reactivated by selecting a new measuring range (4) or a new operating mode (5). See also section 5.2 "Buzzer".

During an overload cut-off, a message concerning the operating instructions will be displayed:



Additionally to the overload cut-off, the instrument is protected by a fuse. A defective fuse can be recognised by the fact that no values are displayed for any of the current measuring ranges. In this case, the fuse must be replaced (see also section 10).

6 FIRMWARE UPDATE

A USB port for service purposes is located under the cover (13). This port can be used to load device updates provided by PHYWE.

6.1 Removing the cover



Attenti-

Ensure that the instrument is switched off and that no measuring leads are connected.

Position the instrument on a level surface with its front panel facing down. On the right side of the back panel's cover, there is a screw. (see Fig. 5)



Fig. 5: Screw holding the cover

Remove the screw with a screwdriver by turning it anti-clockwise. Then, remove the cover.

7 TECHNICAL DATA

7.1 Measuring ranges

Typical at 25°C, warm-up period of 15 min, in battery mode

Selector / unit	Measuring range	Input resistance (approx.)	Accuracy		
			— (DC)	~ (AC)*	
V	1; 3; 10; 12; 30; 100; 300	10 MΩ	±1,5 % EV	±2 % EV	and for: 1 ... 10kHz 2% MV*** 10 ... 100kHz 30% MV
mA****	1	350 mΩ	±1,5 % EV	±3 % EV	
	3; 10; 30; 100; 300			±2 % EV	
A	1; 3; 10	45 mΩ	±1,5 % EV	±2 % EV	

* From 15% of the measuring range end value. To reduce the error, change to smaller measuring range, reference condition 50 Hz

** EV – from measuring range end value

*** MV – from measuring value

**** Depending on the accuracy of the measuring range and the ambient temperature, the device may require a warm-up time of more than 15 min.

6.2 Firmware update

After the cover has been removed (see section 6.1 "Removing the cover"), the USB port (14) for the connection of a PC is accessible.

The firmware and the scales can be updated with the aid of the PHYWE ADM 3 Update Program.

Available updates can be found at:

www.phywe.de >> Downloads >> Softwaredownload

Please check whether the update is more recent than the firmware version of your instrument. Otherwise, an update is not necessary. The version number of your instrument will be displayed briefly when the instrument is switched on. To perform an update, open the program and follow the instructions on the screen of the PC.

After the firmware update, remove the USB cable, place the cover on the back panel, again, and fix it with the screw.



Attention!

Do not use the instrument unless the cover is closed.

7.2 General

Display	TFT display
Digits high	Analogous: 15 mm Digital: 27 mm
Measuring rate	1000 measurements / s
Position of use	upright
Measuring category	CAT II, 300 V max.
Electrical safety	DIN EN 61010-1
Rectifier principle (⎓)	Root-mean-square value (true RMS-to-DC converter)
Frequency bandwidth (⎓)	2 Hz... 100 kHz
Signal shape	any
Waveshape error at a crest factor ≤4	≤0,5 % of the indicated value
Setting time ~	<4 s
Maximum load capacity of the measuring ranges	
Voltage	300 V
Current	10 A
Power supply	2x lead gel battery 6 V, 3,4 Ah
Battery life	approx. 10 hours
Supply voltage	12 V DC, 2,5 A max.
Power-on time	4 s
Dimensions (mm)	380 x 314 x 120
Weight	4,4 kg
<u>Power supply</u>	
Supply voltage	Input 100...240 V AC, 50...60 Hz Output 12 V DC, 2,5 A
Other data	see the operating instructions to article 11262-99

7.3 Ambient conditions

Area of use	Indoor
Operating temperature	5 °C...40 °C
Humidity	max. 80 % rel. hum. up to 31 °C, linearly decreasing 50 % rel. hum. at 40 °C
Operating altitude	2000 m msl max.

8 SCOPE OF SUPPLY

The article is supplied with the following components:

- PHYWE Demo multimeter ADM 3 light
- Power supply 12 VDC/2.5 A 11262-99

Only use original PHYWE spare parts. For replacement, turn to PHYWE or an authorized dealer:

- Overload fuse (1 piece)
FF 10 A, 500 V SP-1384001

9 ACCESSORIES

The following accessories are available:

Measuring leads

- Safety connecting cables, e.g. 07337-88

10 REPLACING THE OVERLOAD SAFETY DEVICE



Attention!

Ensure that the instrument is switched off and that no measuring leads are connected.

Prior to replacing the fuse, ensure that the cause for the blown fuse is eliminated. Never bypass or repair fuses. Open the cover of the fuse holder (12) by turning it anti-clockwise with a screwdriver.

Then, the overload safety device can be replaced and the cover can be closed.

Fuse type: Type G, 10.3 mm x 38 mm
FF 10 A, 500 V, breaking capacity: >10 kA
Application category: gR



Attention!

Do not use the instrument unless the fuse holder is closed!

Should this fuse blow when the instrument is switched on, never replace it with a more resistant fuse! A defect is indicated and the instrument must be returned to the PHYWE service department for repair.

11 WARRANTY

We give a warranty of 24 months for units that we have supplied inside the EU, and a warranty of 12 months outside the EU. The following is excluded from the warranty: damage that is due to non-compliance with the operating instructions, improper use or natural wear.

The manufacturer can only be held liable for the function and safety-relevant properties of the unit if the maintenance, service and modifications of the unit are performed by the manufacturer or by an institution that is expressly authorised by the manufacturer.

12 DISPOSAL

The packaging mainly consists of environmentally-friendly materials that should be returned to the local recycling stations.



Do not dispose of this product with normal household waste. If this unit needs to be disposed of, please return it to the address that is stated below for proper disposal.

PHYWE Systeme GmbH & Co. KG
Customer Service
Robert-Bosch-Breite 10
37079 Göttingen
Germany
Telephone +49 (0) 551 604-274
Fax +49 (0) 551 604-246

13 CLEANING

Only use a soft, dry cloth for cleaning the instrument. The ingress of moisture into the openings of the instrument must be absolutely prevented.

14 EXPLANATION OF THE SYMBOLS



Attention!

Potentially harmful situation
(property damage)
General hazard area
Compliance with the operating instructions is required



Important information about the device
Compliance with the operating instructions is required



Direct current and alternating current



Continuous double insulation

TRMS

True root mean square

CAT II

Measuring category II
(see chapter 2)