# Digital Lock-In Amplifiers

SR850 — DSP lock-in amplifier with graphical display



SR850 DSP Lock-In Amplifier



- · 1 mHz to 102.4 kHz frequency range
- · >100 dB dynamic reserve
- · 0.001 degree phase resolution
- · Time constants from 10 µs to 30 ks (up to 24 dB/oct rolloff)
- · Auto-gain, -phase, -reserve and -offset
- · Data logging (up to 65k samples)
- · Smoothing, curve fitting & statistics
- · GPIB & RS-232 interfaces

to 512 Hz, you are able to see exactly how your data changes in time—not just what the current output value is. After the data has been acquired, the SR850 offers a variety of data reduction options, such as Savitsky-Golay smoothing, curvefitting and statistical analysis. A built-in 3.5" disk drive, along with standard RS-232 and GPIB interfaces, makes it easy to transfer data to your computer.



Large numeric readout with bar graph

The SR850 is a digital lock-in amplifier based on an innovative DSP (Digital Signal Processing) architecture. The SR850 boasts a number of significant performance advantages over traditional lock-in amplifiers—higher dynamic reserve, lower drift, lower distortion, and dramatically higher phase resolution. In addition, the CRT display and 65,536 point memory make it possible to display and process data in a variety of formats unavailable with conventional lock-ins. **Digital Precision** 

At the input of the SR850 is a precision 18-bit A/D converter which digitizes the input signal at 256 kHz. The A/D converter, together with a high-speed DSP chip, replace the analog demodulator (mixer), low-pass filters and DC amplifiers found in conventional lock-ins. Instead of using analog components, the SR850 is implemented by a series of precise mathematical calculations which eliminate the drift, offset, non-linearity and aging inherent in analog components. The same DSP chip digitally synthesizes the reference oscillator, providing a source with less than — 80 dBc distortion, 100 mHz frequency resolution, and 2 mV of amplitude resolution. Digital Flexibility

The SR850 has a 7" CRT display which supports a large selection of display options. Data can be viewed numerically or graphically in bar graph, polar plot and strip chart formats. With 65,536 points of memory and data acquisition rates up

## **Input Channel**

The SR850 has a differential input with 6 nV/ $\sqrt{\text{Hz}}$  input noise. The input impedance is 10 M $\Omega$ , and minimum full-scale input voltage sensitivity is 2 nV. The input can also be configured for current measurements with selectable current gains of  $10^6$  and  $10^8$  V/A. A line filter (50 Hz or 60 Hz) and a  $2\times$  line filter (100 Hz or 120 Hz) are provided to eliminate line related interference. However, unlike conventional lock-in amplifiers, no tracking band-pass filter is needed at the input of the SR850. This filter is used by conventional lock-ins to increase dynamic reserve. Unfortunately, band-pass filters also introduce noise, amplitude and phase error, and drift. The DSP based design of the SR850 has such inherently large dynamic reserve that no tracking band-pass filter is needed.

## **Reference Channel**

The reference source for the SR850 can be an externally applied sine wave or square wave, or its own digitally synthesized reference source. Because the internal reference source is synthesized from the same digital signal that is used to multiply the input, there is virtually no reference phase noise when using the internal reference. The internal reference can operate at a fixed frequency or can be swept linearly or logarithmically over the entire operating range of 1 mHz to 102.4 kHz. Harmonic detection can be performed at any integer harmonic of the reference frequency—not just the first few harmonics.

The DSP approach also offers considerable advantages when working with an external reference. The time to acquire an external reference is only 2 cycles + 5 ms (or 40 ms, whichever is greater)—about ten times faster than conventional lock-ins.

Because the SR850 uses a digital phase-shifting technique rather than analog phase-shifters, the reference phase can be adjusted with one millidegree resolution. In addition, the X and



# Y outputs are orthogonal to within one millidegree. **Outputs** and Time Constants


The output time constants on the SR850 are implemented digitally. Low-pass-filter rolloffs of 6, 12, 18 and 24 dB/octave are available, with time constants ranging from 10  $\mu$ s to 30 ks. Below 200 Hz, the SR850 can perform synchronous filtering. Synchronous filters notch out multiples of the reference frequency—an especially useful feature at low frequencies where the proximity of the 2f component would otherwise require a long time constant for effective filtering. The SR850 makes working at low frequencies a far less time consuming task. **High Dynamic Reserve** 

The dynamic reserve of a lock-in amplifier at a given fullscale input voltage is the ratio (in dB) of the largest interfering signal to the full-scale input voltage. The largest interfering signal is defined as the amplitude of the largest signal at any frequency that can be applied to the input before the lock-in cannot measure a signal with its specified accuracy.

The SR850 has the highest dynamic reserve (>100 dB) of any lock-in available. In conventional lock-in amplifiers, dynamic reserve is increased at the expense of stability. Because of the digital nature of the filtering and gain process in the SR850, the ultra-high dynamic reserve is obtained without any sacrifice in stability or accuracy. In addition, the SR850's high dynamic reserve is obtained without the use of analog band-pass filters, eliminating the noise and error that such filters introduce.

#### **Traces and Displays**

Data acquired by the SR850 is stored in up to four userdefined traces. Each trace can be configured as (A  $\times$  B) / C, where A, B and C are selected from X, Y, R,  $\theta$ , noise, frequency or any of the four rear-panel auxiliary inputs. Common operations, such as ratioing, can be performed in real time by

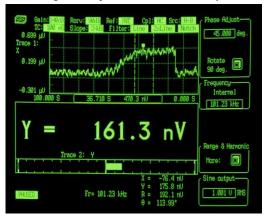

Graphical, numerical, and bar-graph display

defining an appropriate trace. Trace values can be displayed as a bar graph with an associated large numerical display, or as a strip chart showing the trace values as a function of time. Additionally, you can display polar plots showing the phasor formed by the in-phase and quadrature components of the



Polar plot display

signal. All displays can be easily scaled from the front panel or over the computer interfaces, and an auto-scale feature is available to quickly optimize the display. The screen can be configured as a single large display, or as two horizontallysplit displays.

#### **Convenient Auto Measurements**

Common measurement parameters are available as singlekey "auto" functions. The gain, phase, dynamic reserve and display scaling can all be set with a single key press. For many measurements, the instrument can be completely configured simply by using the auto functions.

# Auxiliary A/Ds and D/As

Four rear-panel A/D inputs allow you to measure external signals with millivolt resolution. The measured values can be incorporated into one of the SR850's trace definitions, or can be displayed on the front panel, or read via either computer interface. Four D/A outputs can provide either fixed output voltages or a voltage level which scans synchronously with

the SR850's frequency scans. Both the A/D inputs and the D/A outputs have a  $\pm 10\ V$  range.

# **Analysis Features**

The SR850's performance doesn't stop once data has been acquired—a full set of data processing features is also included. Multiple-range Savitsky-Golay smoothing can be applied to any of the trace arrays, and statistical information (mean, variance, sum) can be calculated for a selected trace region. A curve fitting routine calculates best fits to lines, exponential curves, and Gaussian curves for any portion of your data. And a trace "calculator" lets you perform a variety of simple arithmetic and trigonometric operations on trace data. Interfaces and Hardcopies



The SR850 comes standard with RS-232 and GPIB interfaces. All instrument functions can be queried and controlled via the interfaces. For convenient debugging, characters received and sent via the interfaces can be viewed on the front panel. Several hardcopy options are available on the SR850. Screens can be dumped to a dot-matrix or LaserJet compatible printer through the standard Centronics printer interface. Displays can also be plotted on any HP-GL compatible plotter via GPIB or RS-232.

Ordering I nformation	
SR850	DSP dual phase lock-in amplifier
SR550	Voltage preamplifier (100 M $\Omega$ , 3.6 nV/ $\sqrt{\text{Hz}}$ )
SR552	Voltage preamplifier (100 $k\Omega$ , 1.4 $nV/\sqrt{Hz}$ )
SR554	Transformer preamplifier (0.091 $nV/\sqrt{Hz}$ )
SR555	Current preamplifier
SR556	Current preamplifier
SR540	Optical chopper



SR850 rear panel



# **Signal Channel**

Voltage inputs Single-ended or differential

Sensitivity 2 nV to 1 V 106 or 108 V/A Current input

Input impedance

Voltage input  $10 \text{ M}\Omega + 25 \text{ pF}$ , AC or DC coupled

Current input 1 k $\Omega$  to virtual ground Gain accuracy  $\pm 1 \% (\pm 0.2 \% \text{ typ.})$ Noise (typ.) 6 nV/√Hz at 1 kHz

> $0.13 \text{ pA/}\sqrt{\text{Hz}}$  at 1 kHz ( $10^6 \text{ V/A}$ )  $0.013 \text{ pA}/\sqrt{\text{Hz}}$  at 100 Hz (10<sup>8</sup> V/A)

Line filters 50/60 Hz and 100/120 Hz (Q = 5) **CMRR** 100 dB at 10 kHz, decreasing by

6 dB/oct above 10 kHz

Dynamic reserve >100 dB (without prefilters)

## **Reference Channel**

Frequency range  $0.001~\mathrm{Hz}$  to  $102.4~\mathrm{kHz}$ Reference input TTL or sine (400 mVpp min.)

Input impedance  $1 \text{ M}\Omega$ , 25 pFPhase resolution  $0.001^{\circ}$ Absolute phase error <1° Relative phase error <0.001°  $90^{\circ} \pm 0.001^{\circ}$ Orthogonality Phase

noise

Int. reference <0.0001° rms at 1 kHz

Ext. reference 0.005° rms at 1 kHz, 100 ms, 12 dB/oct

Phase drift <0.01°/°C below 10 kHz,

 $<0.1^{\circ}/^{\circ}$ C, 10 kHz to 100 kHz 2F, 3F, ... nF to 102.4 kHz

Harmonic detection Acquisition time (2 cycles + 5 ms) or 40 ms,whichever is greater

# **Demodulator**

Stability

Digital outputs no drift

Analog outputs <5 ppm/°C for all dynamic reserves

Harmonic rejection -90 dB

Offset/Expand  $\pm 100$  % offset, expand up to 256× Time constants 10 µs to 30 ks (6, 12, 18, 24 dB/oct

rolloff)

Sync. filtering available below 200 Hz

# **Internal Oscillator**

Range 1 mHz to 102.4 kHz Accuracy  $25 \text{ ppm} + 30 \mu Hz$ Resolution 0.01 % or 0.1 mHz (whichever is greater) Distortion

-80 dBc (f < 10 kHz)

-70 dBc (f > 10 kHz) at 1 Vrms

Amplitude 0.004 Vrms to 5 Vrms into 10  $k\Omega$ 

(2 mV resolution)

Output impedance  $50 \Omega$ Amplitude accuracy 1 % Amplitude stability 50 ppm/°C

Sine and TTL (both can be phaselocked to an external reference)

Sweeps Linear and log Stanford Research Systems

**Inputs and Outputs** 



Interfaces IEEE-488.2, RS-232 and Centronics

interfaces standard. All instrument functions can be controlled and read

though the interfaces.

 $X, Y ext{ outputs}$   $\pm 10 V, updated at 256 ksamples/s$  CH1 output  $\pm 10 V ext{ output of } X, R ext{ or Trace 1 to 4}$  CH2 output  $\pm 10 V ext{ output of } Y, \theta ext{ or Trace 1 to 4}$  4 BNC inputs, 1 mV res.,  $\pm 10 V ext{ 4 BNC outputs}$  4 BNC outputs, 1 mV resolution,

±10 V (fixed or swept amplitude) Internal oscillator analog output Internal oscillator TTL output

TTL out Internal oscillator TTL output
Trigger In TTL signal starts internal oscillator
sweep or triggers instrument

data taking (rates to 512 Hz).

Remote pre-amp Provides power to the optional

SR55X preamps

# **Displays**

Sine out

Screen format Single or dual display

are defined as  $A\times B$  / C or  $A\times B$  /

 $C^2$ 

where A, B, C are selected from X,

Y,

R,  $\theta$ , X-noise, Y-noise, R-noise,

Aux 1 to 4 or frequency.

polar plot and strip chart

Data buffer 64k data points. Buffer is configured

as a single trace with 64k points, two traces with 32k points each, or four traces with 16k points each. 0.0625 to 512 Hz, external to 512 Hz

# **Analysis Functions**

Sample rate

Smoothing 5, 9, 17, 21, 25 pt. (Savitsky-Golay)
Curve fitting Linear, exponential or Gaussian
Calculator Arithmetic, trigonometric and

logarithmic calculations
Mean and standard deviation

## General

Statistics

Hardcopy Screen dumps to dot-matrix or

LaserJet printers. Plots to HP-GL

plotters (RS-232 or GPIB).

Disk drive USB drive. Storage of data and

instrument setups (binary or ASCII). Screens can be saved as PCX files.

Power 60 W, 100/120/220/240 VAC,

50/60 Hz

Dimensions  $17" \times 6.25" \times 19.5"$  (WHD)

Weight 40 lbs.

Warranty One year parts and labor on defects

in materials and workmanship



