

Dino-Lite Premier 5018 Series

Dino-Lite
Digital Microscope

NOTICE

- It is recommended to use Dino-Lite in a well-ventilated environment.
- Be sure to use the supplied power adaptor connecting to Dino-Lite for preventing any damage to the device.


Package Contents

- Dino-Lite Premier
- DVI cable
- Power adaptor (5V)
- Carrying pouch


Features

- Direct digital DVI interface for LCD or HDTV monitors.
- Sharp and vibrant HD 720p resolution.
- Dual purpose MicroTouch to freeze the frame or switch LEDs.
- Adjustable magnification varied by model:
 - AM5018 series: 20x-50x, 200x
 - AD5018 series: 20x~200x
 - AM5018 L / AD5018 L series: 20x~90x

System Configuration

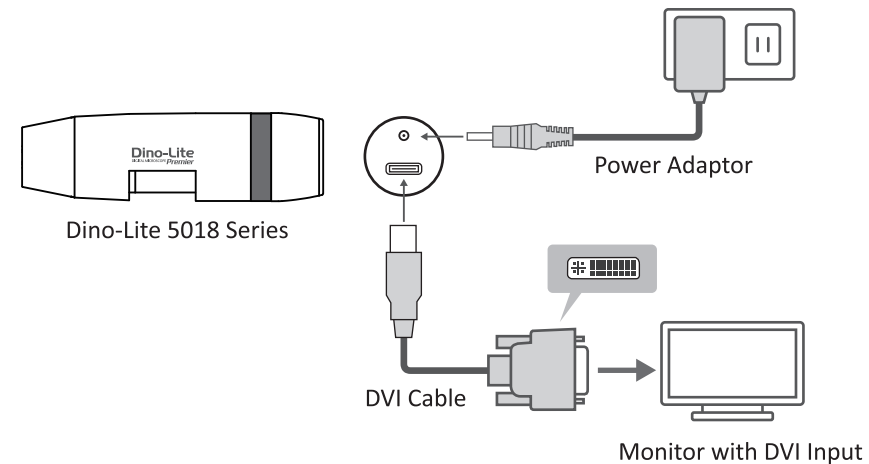

How to Use

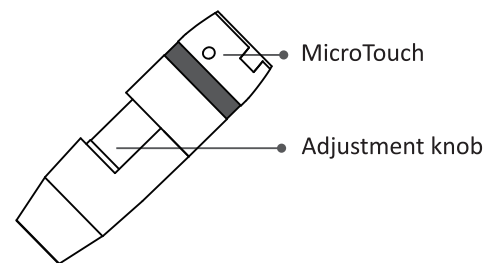

Location of Dino-Lite Controls

- Use the *adjustment knob* to obtain a focus or magnification with the Dino-Lite.
- Tap the *MicroTouch* to freeze / unfreeze frame.
- Tap and hold the *MicroTouch* for 2 seconds to switch the LEDs.

For more information, please visit: www.dino-lite.com

Made in Taiwan



P/N:2410181