WHM

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Contents

1 Introduction	8
2 Safety Instructions	8
2.1 Description of symbols	8
2.2 Instructions	8
3 Product description	10
3.1 Key features	10
3.2 General survey	10
3.3 Lamps and control elements	10
3.4 Restrictions	10
4 Package content	11
5 Positioning the device and commissioning	11
6 Job description	11
7 Maintenance and cleaning	11
8 Technical data	12
9 Certificate	13

1 Introduction

Thank you for purchasing our **WHM Multi Heating Mantle with analog Controller**. You did a good choice buying a **Wisdom** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Due to its functions the Heating Mantle is not protected against incoming water. Only a professional handling can avoid hazardous accidents.

- Only persons with sufficient safety knowledge are allowed to use this device.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Do not touch the device if any liquid is flowed into the device (e.g. broken glass vessel) and firstly switch the power supply off. After that please pull out the power plug.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Do not put heat-sensitive items at the device.
- Only use adequate sized vessels and avoid an over boil of the liquids.
- Do not clean the scorching device.

- Use this device only if you read the whole safety instructions.
- This device should only used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.
- Please make sure that you use only vessels with heat resistance glass (look at 3.4).
- Do not heat up the heating mantle without any vessel to avoid an overheating.
- Do not heat the up the device more than it is necessary - for lower energy consumption and your own safety.

- Make sure that the device can not slip away.
- To avoid heat congestions or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in areas with potentially explosive atmosphere.


- Repairs and amendments should only be made by **witeg** authorized people. Contraventions exclude the warranty.

3 Product description

3.1 Key features

The **Wisdom WHM Multi Heating Mantle** is designed for chemical, biological, medical and pharmaceutical heating and extraction applications up to 450 °C. We offer devices with 3 and 6 heating places for round flasks.

3.2 General survey

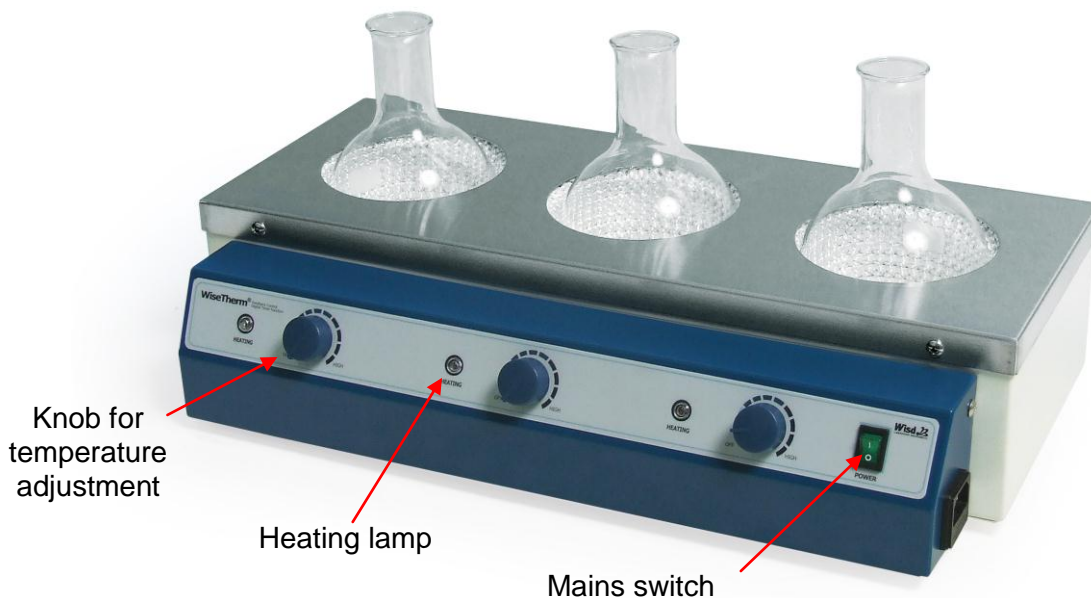Knob for temperature adjustment

Heating lamp

Mains switch

3.3 Lamps and control elements

Knob for temperature

- Adjustment of temperature.
- Until left stop: heater is off.
- Until right stop: heater is on.

Heating lamp

- Indicates heating process.

Main switch

- On / Off switch.

3.4 Restrictions

Please only use glass vessels that are made for heating applications. A minimum heat resistance like borosilicate glass 3.3 has is needed.

4 Package content

quantity	denomination
1	WHM Multi Heating Mantle with analog Controller
1	operating manual

5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your costumer service.
4. Remove perhaps soiling and package material.
5. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
6. Make sure that you follow all safety instructions.
7. Connect the device with the power outlet.
8. Switch the device on and start the process (look at job description).

6 Job description

Set the temperature by turning the rotary knob. Turn it until the left stop to switch out the heating. The heating lamp shines if the heating is activated.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts and accessories.
- Make sure that no liquid gets into the device.

8 Technical data

Multi Heating Mantle							
Temperature		Max. 450 °C					
Material Heating Area		Glass cloth					
Material Housing		Powder coated Aluminium					
Safety class		IP10					
Article number DH.WHM12...		Heating places	Round flask size (ml)	Flask Ø (mm)	Power (W)	Inside depth (mm)	Housing size (mm) W x D x H
230 V 50 Hz	110 V 60 Hz						
391	392	3	500	102-105	720	65	550 x 270 x 130
393	394	3	1.000	130-133	1.050	75	690 x 340 x 170
291	292	6	230/300	83-86	840	50	800 x 240 x 110
293	294	6	250/300	83-86	840	50	800 x 240 x 110
295	296	6	500	102-105	1440	65	950 x 270 x 130

9 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	WHM Multi Heating Mantle with analog Controller
Typ	
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
 D-97877 Wertheim
 TEL: +49-9342-9301-0
 FAX: +49-9342-9301-77
 Email: info@witeg.de
 www.witeg.de