WIS-ML

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Multi-Stack Shaking Incubator

Operating Manual

gilt für / written for
WIS-ML02/ -ML04

witeg Labortechnik GmbH

(English manual at page 13)

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Contents

| | |
|---|-----------|
| 1 Introduction | 14 |
| 2 Safety Instructions | 14 |
| 2.1 Description of symbols..... | 14 |
| 2.2 Instructions | 14 |
| 3 Product description | 16 |
| 3.1 Key features..... | 16 |
| 3.2 Operating surface..... | 17 |
| 3.3 Term description | 17 |
| 4 Accessories | 18 |
| 5 Positioning the device and commissioning | 18 |
| 6 Job description..... | 19 |
| 6.1 Temperature Settings..... | 19 |
| 6.1.2 Set turn-on delay | 19 |
| 6.1.3 Set temperature | 19 |
| 6.1.4 Justify | 20 |
| 6.1.5 Key lock | 20 |
| 6.1.6 Storage function | 20 |
| 6.2 Shaking process | 21 |
| 6.2.1 Menu structure | 21 |
| 6.2.2 Set speed and delay timer | 21 |
| 7 Maintenance and cleaning..... | 22 |
| 8 Troubleshooting Guide..... | 22 |
| 9 Technical data..... | 23 |
| 10 Certificate..... | 24 |

1 Introduction

Thank you for purchasing our **WIS-ML Multi-Stack Shaking Incubator**. You did a good choice buying a **Wisdom** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**Risk of injury**

Avoid injuries of your hands by following these instructions.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Ensure that the power supply cord is undamaged before you connect it with the power source.
- Do not use power bars to extend the power supply cord.
- Use the device only in electric circuits with ground fault circuit interrupter (RCD). The same applies to devices with integrated fault circuit switch.
- Only use extension cables with sufficient cable cross section.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- If easy flammable, electro conductive or other objects reached the inner of the devices housing switch out the device immediately and disconnect it the electrical power source.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- It is important that the device stands up on a horizontal, stabile and firm subsurface to avoid vibrations.
- Make sure that the device can not slip away.
- The device is easy moveable with its rollers and can be easily fixed at its working place. Therefore turn the red wheels at the rollers to pull down the stands.
- Make sure that the device is standing safely and equally at all its stands.
- To avoid heat congestion or fires make sure that there is enough space for air circulations. Therefore is a minimum distance of 20 cm needed between ventilation openings and next objects.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in area with potentially explosive atmosphere.

- Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.
- Make sure, that the safety switches are adjusted up to 30 °C over maximum and minimum temperature. Otherwise the set temperature may be not reached.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- Wear corresponding protective work clothing during the usage.
- Switch the device off before you disconnect it from the electrical power source.
- Pull out the power cable if you didn't want to use the device for a longer time.
- You have to keep the device clean if you want to get a longer life of it.


- Do not put your hands into the device as long as the shaking mechanism works.
- Pull out the power supply cord before you clean the device.

3 Product description

3.1 Key features

- Long life by simply and solid construction.
- Shaking speed up to 250 rpm and separately adjustable for each stack
- Turn-off and turn-on delay timer with signal sound
- Temperature range from 10 °C up to 60 °C
- Door window
- Quiet running
- Automatic open door stop
- Easy to justify
- Key lock against inadvertent changing of set temperature and time values
- Illuminated LCD
- Storage function for adjusted time and temperature values.
- Universal platform for flask holders included
- Shaking rack available (optional accessories).

3.2 Operating surface

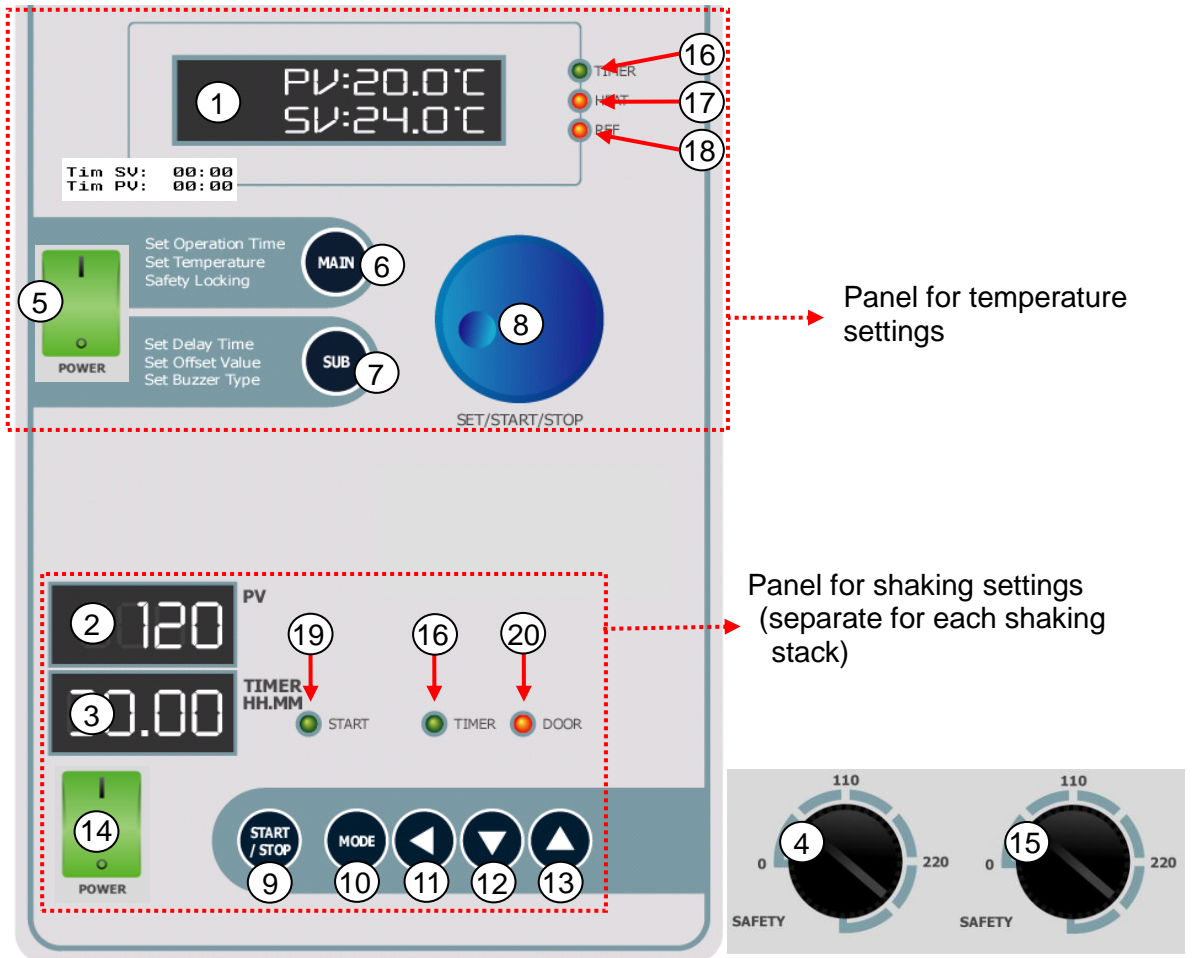3.3 Term description

- 1: LCD (Temperature):** Illuminated LCD for temperature controller
 - Upper line shows set value (SV) and lower line present value (PV).
- 2: PV Display (Shaking):** display for actual value, PV = present value
- 3: SV Display (Shaking):** display for Timer and set value, SV = set value
- 4: Over temperature safety switch**
 - The safety switch turns the heating off if your set temperature is overshoot.
 - Set the maximal temperature rough 20 – 30 °C over the desired maximal safety temperature, otherwise the device maybe not reaches your set value.
- 5: Main switch for temperature controller**
 - On / off switch for temperature controller.
- 6: MAIN button**
 - For changing between the different menu items.
- 7: SUB button**
 - For reaching the sub menu Items.
- 8: Jog-Shuttle knob**
 - Change SV by turning.
 - Push the knob for start and stop or probate set values.
 - Hold id pushed for two seconds to save set values.
- 9: Start/Stop button (Shaking)**

10: MODE button (Shaking)

- Change between speed, Timer and keylock menu items.

11: Button left arrow (shaking)

12: Button right arrow (shaking)

13: Button up arrow (shaking)

14: Main switch for shaking controller

- On / off switch for shaking controller.

15: Over temperature safety switch

- The safety switch turns the heating off if your set temperature is undershoot.
- Set the maximal temperature rough 20 – 30 °C over the desired minimal safety temperature, otherwise the device maybe not reaches your set value.

16: Timer lamp

- Is shining when turn-off delay timer runs.

17: Heating lamp

- Is shining when heater works.

18: Refrigerating lamp

- Is shining when refrigerator works.

19: Shaking lamp

- Is shining when shaking mechanism works.

20: Door lamp

- Shines when door is open.

4 Accessories


| Quantity | Denomination |
|----------|---|
| 1 | WIS-ML Multi-Stack Shaking Incubator |
| 1 | Operating manual |

5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol: 
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your costumer service.
4. Make sure that you follow all safety instructions.
5. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
6. Ensure that the device was not tilted more than 20 ° within the last 24 hours. If you can not be sure so please let it stand upright for 24 for hours before first usage. Otherwise the cooling unit may be damaged.
7. Turn the red wheels to lower the stands. That will fix the device at its working ground.
8. Set the safety switches (3.3).
9. Connect the device with the power outlet.
10. Switch the ground fault circuit interrupter on.
11. Switch the device on.
12. Set your values and start the process (look at job description).

6 Job description

6.1 Temperature Settings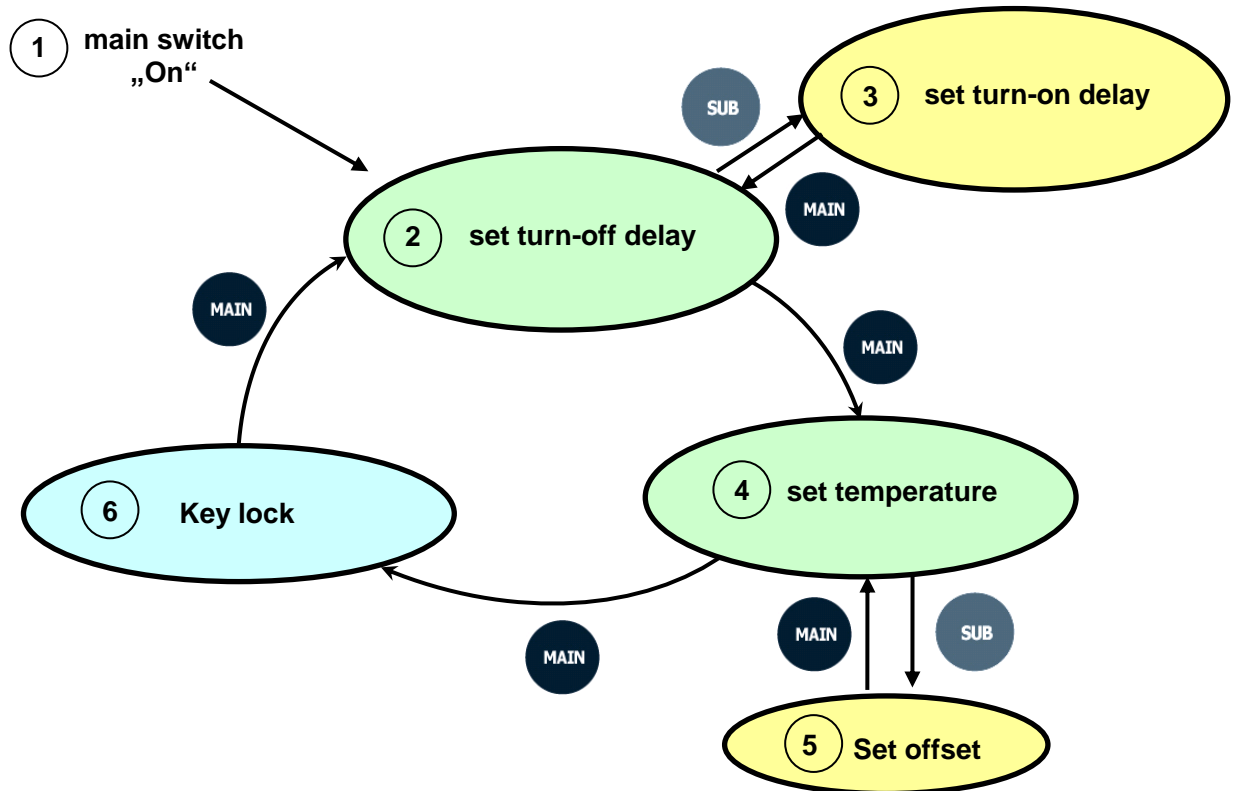

6.1.1 Set turn-off delay

The turn-off delay timer serves as automatically stop function. It turns off the heating and pump automatically.

- First menu item after turning on is “set turn-off delay”. The timer lamp is blinking.
- Turn the Jog-Shuttle to set the desired turn-off time.
- Start and stop timer by pushing the Jog-Shuttle knob.
- Timer must be stopped to enable a renewed changing of timer value.
- In the upper display line is the SV (set value) and in the lower line the remaining time to turn-off (PV) shown.
- Stop timer for endurance run.

6.1.2 Set turn-on delay

The turn-on delay timer serves as automatically start function. It turns on the heating and pump. It is possible to work with turn-off timer in addition.

- Have a look to the menu overview (6.1) to open the relevant menu item.
- Turn the Jog-Shuttle to set the desired turn-on time.
- Start and stop timer by pushing the Jog-Shuttle knob.
- Timer must be stopped to enable a renewed changing of timer value.
- In the upper display line is the SV (set value) and in the lower line the remaining time to turn-on (PV) shown.
- Stop timer for immediately run.

6.1.3 Set temperature

- Have a look to the menu overview (6.1) to open the relevant menu item.

- Turn the Jog-Shuttle to set the desired temperature.
- Start and stop heater by pushing the Jog-Shuttle knob.
- Heater must be stopped to enable a renewed changing of temperature.
- In the upper display line is the SV (set value) and in the lower line the actual reached temperature (PV) shown.
- Stop timer for immediately run.

6.1.4 Justify

May the device doesn't show the real temperature you can justify it.

- Calculate the offset as follows:

$$\text{Real temperature} - \text{shown Temperature (PV)} = \text{Offset}$$

- Have a look to the menu overview (6.1) to open the relevant menu item.
- Turn the Jog-Shuttle to set the desired offset
- Push the main button to confirm the offset. Thereby you reach the menu item "set temperature".

6.1.5 Key lock

- To avoid accidental changing of the set values, push the MAIN button in the menu item "set temperature".
- In this menu item can you see the actually temperature in the SV-line and the remaining time of turn-off delay in the PV-line. The function of the Jog-Shuttle knob is deactivated.
- Press the MAIN button once again to leave the key lock. Than you are in the menu item "set turn-off delay".

6.1.6 Storage function


Storage Function is very useful if you want to work with the same parameters often. Then you didn't need to set them in every usage.

- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV for three times.
- Any stored parameters are erasable. Turn the value with the Jug-Shuttle till "0" and then hold the knob pressed for two seconds.
- If you want to overwrite a value it is not necessary to erase the old one. Set the desired value and use the Jog-Shuttle knob for storing.

6.2 Shaking process

6.2.1 Menu structure

All settings for shaking are made with buttons 9 to 13. Display 2 and 3 are for visualisation.

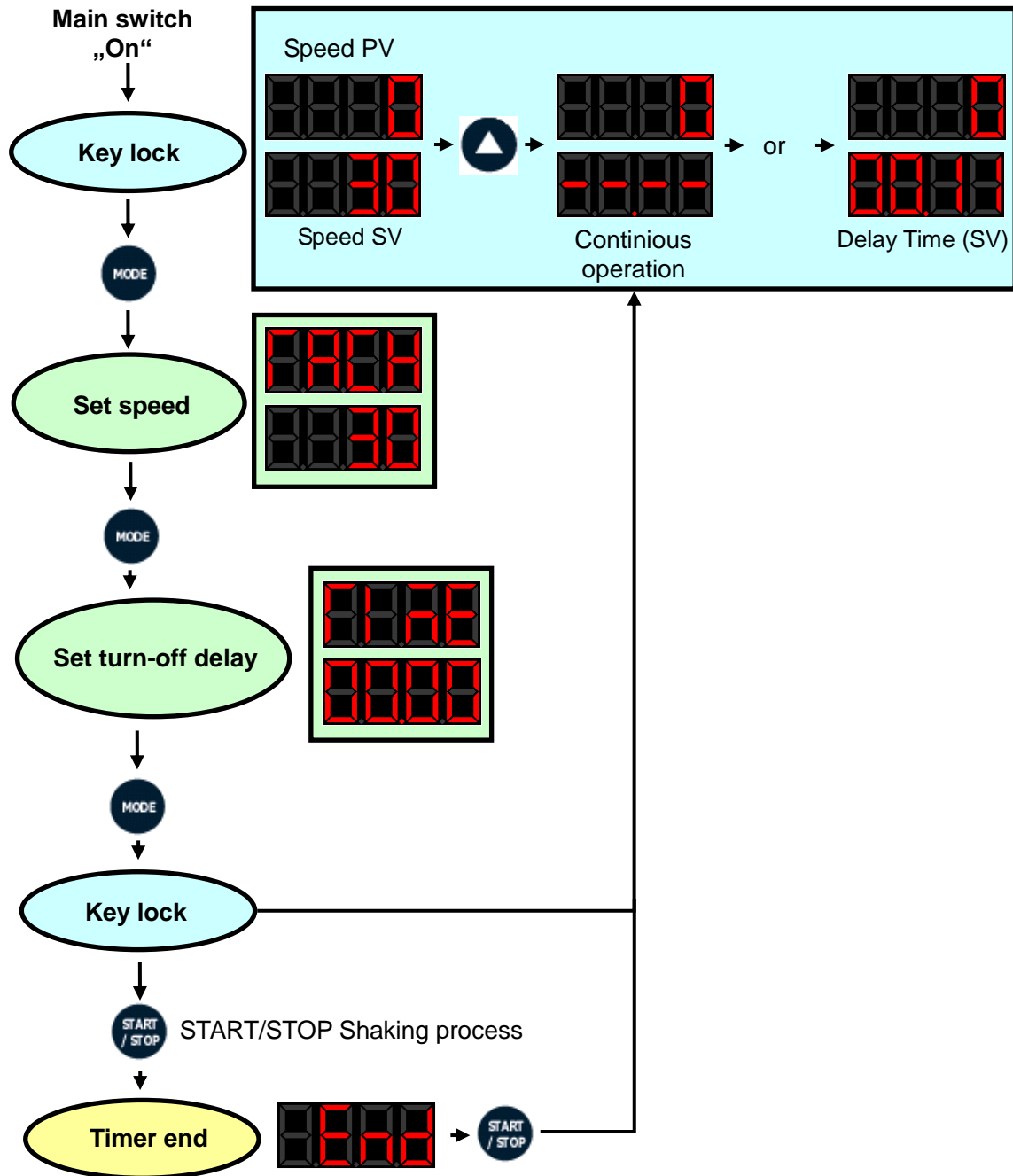6.2.2 Set speed and delay timer

Have a look to the menu structure (6.2.1) to reach the menu items for timer and speed setting. When you enter a menu item the last digit starts blinking. With the buttons 12 and 13 you can change the values of the blinking digits. Use button 11 to choose the other digits.

When you push the MODE button the set value is stored and you change to the next menu item.

The timer menu item has a dot behind the second digit. The digits before the dot are for hour setting and behind the dot for minutes.

To start and stop the shaking process you have to enter the key lock menu item. Push Start/stop button for start and stop.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry.
- Only use original spare parts.
- Check the cleanliness of the fin condenser in regular intervals of 6 month. You can vary this periods depending to the air pollution. The fins have to be as clean that the heat easily can be blown away by the fan. You can use pressure air, brushes or vacuum cleaners. Do not bend the fins!

8 Troubleshooting Guide

| Situation | Confirmations and Solutions |
|---|---|
| Power Failure | <ul style="list-style-type: none"> ✓ Check the supplied voltage is present. ✓ Check the power connection and the power cord for proper connection and integrity. ✓ Check if the MAIN Power Switch is "ON". ✓ Fuse is blown. |
| Unit Not Heating | <ul style="list-style-type: none"> ✓ Check if the MAIN Power Switch is "ON". ✓ Check if the set temperature (SV) confirmed with the Jog-Shuttle knob (by turn the Jog-Shuttle knob cannot change the set temperature – see point 6.1.3). ✓ Check the Safety switch is set higher (min. 10°C-20°C) than the set temperature (SV). |
| Heating, but unit not reaching temperature or temperature is fluctuated | <ul style="list-style-type: none"> ✓ Check the Safety switch is set higher (min. 10°C-20°C) than the set temperature (SV). ✓ Problem in the sensor or main controller. |
| Unit Not Refrigerating | <ul style="list-style-type: none"> ✓ Check if the MAIN Power Switch is "ON". ✓ Check if the set temperature (SV) confirmed with the Jog-Shuttle knob (by turn the Jog-Shuttle knob cannot change the set temperature – see point 6.1.3). ✓ Check that the set temperature (SV) is lower than 35 °C. |
| Refrigerating but it doesn't fall to the SV temperature | <ul style="list-style-type: none"> ✓ Problem with the the cooling circuit. |
| Unit Not Shaking | <ul style="list-style-type: none"> ✓ Check if the MAIN Power Switch is "ON". ✓ Check if the set RPM confirmed with the Start button. |
| ERROR 1 | <ul style="list-style-type: none"> ✓ Temperature Sensor or main controller Error. |
| ERROR 3 | <ul style="list-style-type: none"> ✓ Door open alarm and appears when you open the door. ✓ If Error 3 develops during door closed, check the Limit Switch is pressed properly. |

If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.

Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.

9 Technical data

| | | WIS-ML02 | WIS-ML04 |
|--------------------------------------|---------------------|--|---------------|
| Dimensions (BxTxH) | Chamber (mm) | 700x700x1100 | 700x700x1790 |
| | Capsule (mm) | 1290x850x1445 | 1290x850x2125 |
| Stacks | | 2 | 4 |
| Heating power | | 1.5 kW | 1.8 kW |
| Cooling power | | 370 W | 550 W |
| Temp. | Range | 10 °C up to 60 °C | |
| | Accuracy | ± 0,1 °C | |
| Shaking | Speed | 30 - 250 rpm | |
| | Amplitude | 25 mm orbital motion | |
| Timer | | 99 h 59 min, endurance run | |
| Refrigerant | | CFC-free (R-134A) | |
| Material | Chamber | Stainless steel | |
| | Capsule | Powder coated steel | |
| Door | | Window, magnet packing | |
| Platform | | Universal platform | |
| Optional accessories | | Spring rack, Flask holder for 50 -1000 ml flasks | |
| Quantity of flask holders | FH100 | 62 ea | |
| | FH250 | 36 ea | |
| | FH500 | 25 ea | |
| Safety devices | | Overheating, undercooling and over current protection | |
| Others | | Memory function (Temp., Timer), Key lock (Jog-Shuttle input disabled) Signal sound (Error and Timer) Automatic open door stop | |
| Power supply | | AC 230 V, 1N~, 50/60 Hz | |

10 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the customer.

Please look on your account and complete following table:

| | |
|-----------|---|
| Article | WIS-ML Multi-Stack Shaking Incubator |
| Typ | |
| Serialno. | |
| Date | |

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de