Incubator Oven

Controller SMART-Lab™ Jog-Dial

Anytime - Anywhere connected laboratory



SMART-Lab™ Controller



- full touch screen TFT LCD:
- WiRe™ app service, remote control system
- variable program settings
- self-diagnosis and interactive temperature graph
- automatic data recording and password protection
- · data transfer to PC by USB memory
- digital calibration (offset function)
- min-/max-temperature memory & program function
- push-alarm service to smartphone or tablet PC

more information on 8 - 11

SMART-Lab devices

Incubator

SWIG 70°C, 32 / 50 / 105 / 155 I - page 59

SWIF 70°C, 50/105/155 I - page 61

SWIR 0-60°C, refrigerated - page 65

STH -20/-40°C, 98% rel. hum. - page 72

STH-E -20°C, 95% rel. hum. - page 73

SWGC Illumination, 95% rel. hum. - page 74

Oven

SWON 230°C, 32/50/105/155 I - page **77**

SWOF 250°C, 50/105/155 I - page **79**


SWOV 200°C, 18.6/30/70 I, 750mmHg - page 81

Jog-Dial controller

- ergonomic design for easy and convenient usage
- 2-way Jog-Dial knob provides simple setting of required temperature and time
- · high quality illuminated LCD
- "MAIN" and "SUB" button to access all menus




Incubator Oven WIG 70°C, 32 / 50 / 105 / 155 I on page 58 WON 230°C, 32/50/105/155 I on page 76 WIF 70°C, 50/105/155 I on page 60 WOF 250°C, 50/105/155 I on page 78 WIR 0-60°C, refrigerated on page 64 WOV 200°C, 18.6/30/70 I, 750mmHg on page 80

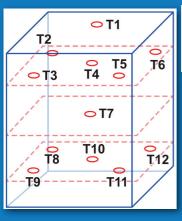


Temperature uniformity successfully tested by **ASTM** standards

All incubators and ovens are thor-

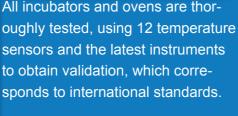

Incubator Oven

General information

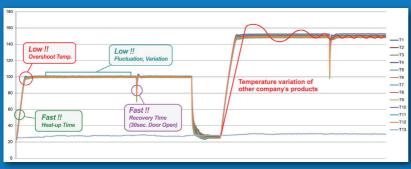

Green product

All incubators and ovens achieve minimum power consumption by using heating elements consuming minimum power adjusted for the capacity and temperature of the chamber.






Incubator air-flow

Gravity-air

The incubators gravity-air flow is established by creating radiant and conductive heat on three sides of the chamber (bottom, left and right side).

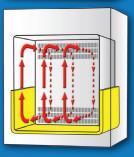


Forced-air

The incubators forced-air flow is established by a three sided heating element in combination with a very effective fan placed in an air-flow room for best temperature uniformity in the whole chamber.

Oven air-flow


Gravity-air

The ovens gravity-air flow is established by heating up air in a pre-heating zone by heating elements on the bottom, left and right side. The pre-heated air will be restrictively provided into the chamber.

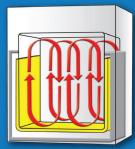

Forced-air

The ovens forced-air flow is established by a very effective fan installed in the rear of the chamber sucking air from the chamber into the pre-heating zone where heating elements on the bottom, left and right side heat up air. The pre-heated air is uniformly and quickly provided into the chamber by a special side and bottom structure.

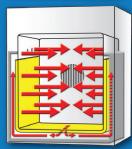

Incubator | growth chamber

Smart-Lab controller, 0-12,000/15,000 Lux, 10°C - 60°C (refrigerated), up to 95% relative humidity









SWGC Illumination, 95% rel. hum.

Ideal for:

- growth of organisms and plants, stability and endurance testing of materials, environmental simulations
- stability product self-life QC testing

Features:

- CFC free refrigerant (R-404A)
- 3-sided light bank system: intensity of light is controlled by 8 levels and displayed on LCD
- inner tempered glass door with silicone packing offers easy observation without door opening
- PE-coated wire shelves included (4 or 8 acc. to model)
- · 3-sided magnetic doors with light banks
- stainless steel interior and powder-coated body
- CE certified and unique serial number for tracing

Safety mechanism:

- push alarm service to your smartphone
- overheat and over-current protection
- sensor error detection and leakage breaker

Controller:

- Smart-Lab™ controller with 7" full touch screen TFT LCD:
- ♦ WiRe™ app service, remote control system
- variable program settings
- ♦ self-diagnosis and interactive temperature graph
- ♦ automatic data recording and password protection
- ♦ data transfer to PC by USB memory
- ♦ digital calibration (offset function)
- ♦ min-/max-temperature memory & program function

Model	SWGC-450	SWGC-1000
Capacity	432 l (1 door)	864 I (2 doors)
Temperature range & variation	10°C - +60°C, ±1°C at 20°C	
Temperature fluctuation	$\pm 0.5^{\circ}$ C at 10°C, $\pm 0.5^{\circ}$ C at 25°C (0% light intensity) / $\pm 1^{\circ}$ C at 10°C, $\pm 1^{\circ}$ C at 25°C (100% light intensity)	
Humidity range, fluctuation & variation	30% - 95% RH, ±1%, ±3% at 60°C	
Humidifier	evaporation by fan with heater	
Illumination lamp & range	14 x FPL 55 W, 0 to 12,000 Lux	28 x FPL 55 W, 0 to 15,000 Lux
Illumination bank	3x light banks: left, right, front	
Temperature & humidity sensor	temperature: PT100, humidity: capacitive sensor	
Heat-up time (up to 37°C)	11 min (0% illumination) / 9 min (100% illumination)	
Cool-down time (down to 10°C)	13 min (0% illumination) / 15 min (100% illumination)	
Recovery time (door open 30sec)	3 min. to 10° C, 1 min. to 37° C (0% illumination); 5 min. to 10° C, 1 min. to 37° C (100% illumination)	
Heating power (temperature & humidity)	1.5 kW	3.0 kW
Compressor	370 W	750 W
Program function	temperature and humidity: 120 patterns, 1.200 segments; light bank: 8 steps	
Shelves	4 x PE-coated steel wire shelves (16 kg / shelf)	8 x PE-coated steel wire shelves (16 kg / shelf)
Internal dimensions (W x D x H)	650 x 700 x 1050 mm	1200 x 790 x 1050 mm
External dimensions (W x D x H)	850 x 1090 x 1965 mm	1200 x 1150 x 1965 mm
Packing size & gross weight	1400 x 1220 x 2210 mm, 435 kg	1600 x 1500 x 2200 mm, 585 kg
Power supply	1 Phase AC 120V, 60 Hz or AC 230V, 50/60 Hz or 3 Phase	
Power consumption	4,8 kW	6,5 kW
Order number 230V*	DH.SWGC00450	DH.SWGC01000
Order number 120V*	DH.SWGC50450	DH.SWGC51000
Order number 3 Phase	DH.SWGC60450	DH.SWGC61000

^{* 230}V and 120V models not available in Europe