
***FP / FHP***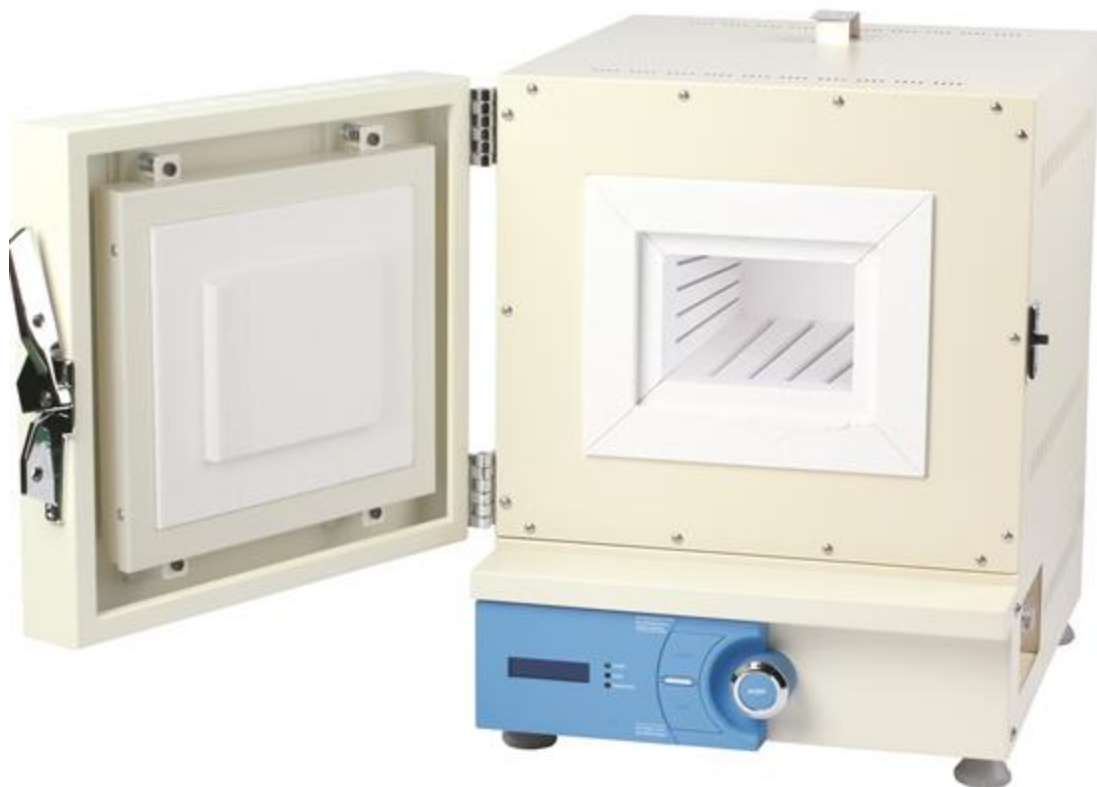

**Programmierbarer Muffelofen**  
**Bedienungsanleitung**  
**Programmable Muffle Furnace**  
**Operating Manual**

gilt für / written for  
FP + FHP -03 / -05 / -12 / -14 / -27 / -63

**witeg Labortechnik GmbH**

(English manual at page 12)

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## 1 Introduction

Thank you for purchasing our **Programmable Muffle Furnace**. You did a good choice buying a **Wisdom** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

**Witeg** reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

## 2 Safety Instructions

### 2.1 Description of Symbols

#### **WARNING**

Information for avoiding injuries or fatal accidents.

#### **ATTENTION**

Safety instruction to avoid a damage of the product.

#### **ATTENTION**

Information about technical parameters.

#### **Environment**

Information about the operating environment.

#### **WARNING of explosions**

Safety instruction to avoid explosions.

#### **ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

### 2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.

- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Because the temperature of the block can reach up to 1,000°C (FP models) or 1.200°C (FHP models), please be careful when opening and closing the Lid.
- Do not put heat-sensitive items at the device.
- Warn all attendees if you leave the device unattended.
- Do not clean the scorching device.

- Use this device only if you read the whole safety instructions.
- This device should only use by persons who are familiarized with safety instructions made for laboratories.
- **Only open the door, if the heating is switched of.**
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.

- It is important that the device stands up on a horizontal, stabile and firm subsurface to avoid vibrations.
- Make sure that the device cannot slip away.
- To avoid heat congestions or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Reparations and amendments should only made by **witeg** authorized people.  
Contraventions exclude the warranty.

- Do not use this device in areas with potentially explosive atmosphere.

### 3 Packing contents

quantum	denomination
1	FP / FHP Muffle Furnace
1	Ceramic Fiber Plate
1	operating manual

## 4 Product description

### 4.1 Key features

The **Programmable Muffle Furnace** is designed for chemical, biological, medical and pharmaceutical heating-applications. The device was developed for simplest usage and a very long life.


Key features are as follows:

- Long life by simply and solid construction.
- 4 side Heating.
- Jog-Dial Knob for easy and fast handling.
- Short heating-up periods.
- Turn-off delay from one minute up to over 99 hours in one minute steps.
- Low weight and high grade insulation.
- Storage functions for set time and temperature values.
- Signal sound after end of turn-off delay and error.
- Key lock against inadvertent changing of set Temperature and timer values.

### 4.2 General survey

### 4.3 Operating surface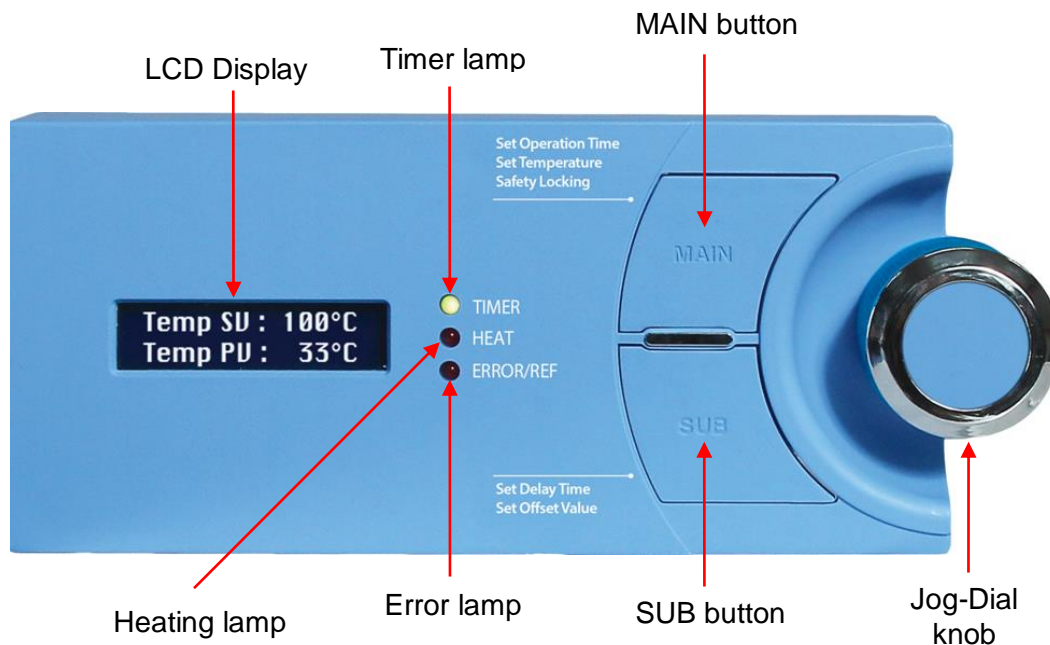

### 4.4 Term description

**SV:** set value (upper line in LCD)

**PV:** present value, measured by the device (lower line in LCD)

#### Main switch

- On / off switch.

#### Timer lamp

- Is blinking in menu items for timer modulation
- Shines continuous in the key lock menu item.

#### Heating lamp

- Shines if the heater is activated.

#### Error lamp

- Is lightning if the door is opened during heating process or another error is appeared.

#### Jog-Dial knob

- Change SV by turning.
- Push the knob for start and stop or probate set values.

#### Safety switch

- The safety switch turns the heating off if your set maximal temperature is overshoot.
- Set the maximal temperature rough 40 – 50 °C over the desired maximal temperature, otherwise the device maybe not reaches your set value.

#### MAIN button

- For changing between the different menu items.

#### SUB button

- For reaching the sub menu Items.

## 5 Positioning the device and commissioning


1. Follow especially the safety instructions marked with following symbol:
2. Remove the package carefully.

3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Remove perhaps soiling and package material.
5. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
6. Make sure that you follow all safety instructions.
7. Connect the device with the power outlet.
8. Switch the device on.
9. Set your values and start the process (look at job description).

## 6 Job description

### 6.1 Menu structure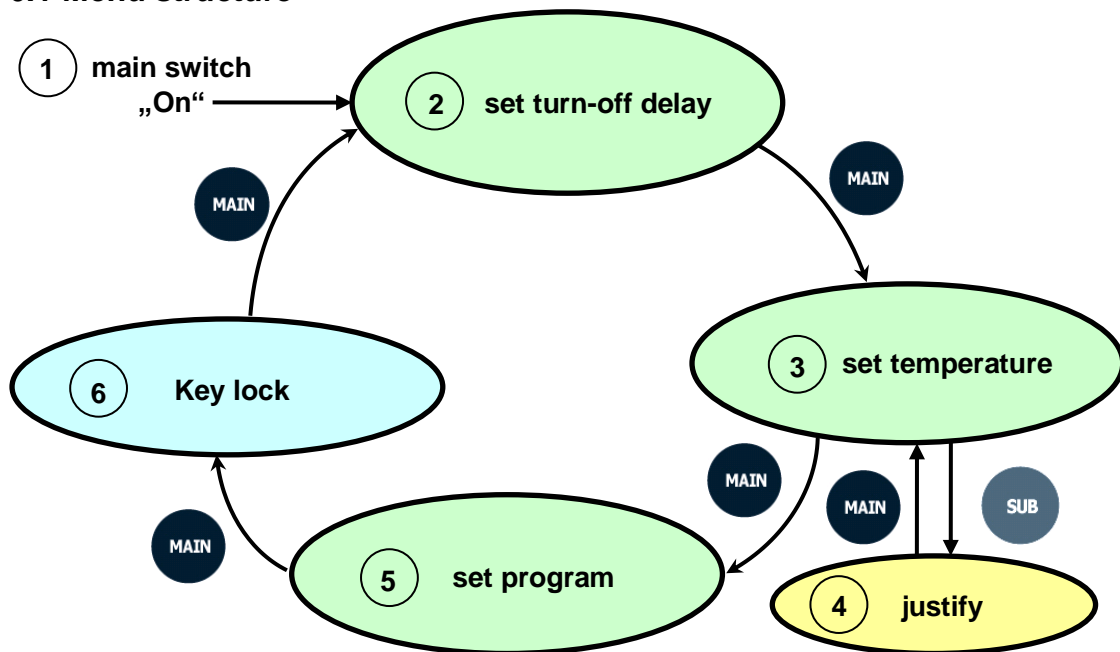

### 6.2 Set turn-off delay

The turn-off delay timer serves as automatically stop function. It turns off the heating after elapsing of the timeframe you set.

- First menu item after switching the device on is the menu “set turn-off delay”. The timer lamp is blinking.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- “00:00” stands for endurance run.
- Turn the Jog-Dial to choose your desired value and probate it by pushing the Jog-Dial.
- If you want to change the confirmed value, you have to push the Jog-Dial once again.

### 6.3 Set temperature

- To reach this menu item you have to push the MAIN button in the menu “set turn-off delay”
- The temperature is adjustable to 1000 °C (F Series) or 1200 °C (FHP Series).
- Turn the Jog-Dial to choose your desired value and probate it by pushing the Jog-Dial. The delay time starts to run by probating the value.
- If you want to change the confirmed value, you have to push the Jog-Dial once again.

- Turn the Dial knob clockwise to increase the temperature and counter-clockwise to reduce it.
- The heating lamp shines, if the heater is activated.

## 6.4 Justify

If the temperature measured by you differs from the temperature the device is showing (PV) you have the possibility to compensate the divergence by the Offset. So you can balance differences of about  $\pm 50^{\circ}\text{C}$ . The device will show the same temperature as the thermometer you used.

The Offset is calculated as follows:

**Temperature measured by you – showed Temperature (PV) = Offset**

- To reach this menu you have to push the SUB button in the menu item “set temperature”.
- You can recognize this menu at the PV showed in the SV-line.
- If you turn the Jog-Dial knob.
- Set the calculated value by turning the Jog-Dial. The Offset will be added to the PV immediately.
- Save and confirm the Offset by pushing the MAIN button, you will also reach the previous menu.

## 6.5 Set Program

- To reach this menu item you have to push the MAIN button in the menu “set temperature”.
- There are 4 Programs (1 ~ 4) available for the setting of each value. And each program allows 10 Steps (0 ~ P) for the setting.

<b>Program no.    1</b> <b>Temp SV : 1000°C</b>
--

- Turn the Jog-Dial-Knob to set the Program No. and push the Jog Dial-Knob for over 1 second to set the STEP.
- Set the STEP of the selected Program No. from the STEP “0” for the values of Set Temperature, Rising Time, and Duration.

<b>STEP :            0</b> <b>Temp SV : 1000 °C</b>
--

- Turn the Jog Dial-Knob to set the value of temperature (Temp SV).
- Push the MAIN button after finishing the setting of temperature to set the value of Rising Time.

<b>STEP :            0</b> <b>Time RV : 01:00</b>
--

- Turn the Jog Dial-Knob to set the value of Rising Time. The Rising Time can be set up to 99 hours 59 minutes; and if the set value of Rising Time is shorter than the output value of the Heater then the program would reach the target temperature with the maximum output value independent of the set Rising Time.


- Push the MAIN button after finishing the setting of the value of Rising Time (Time RV) to set the value of Duration (Time MV).

**STEP : 0**  
**Time MV : 02:00**

- Turn the Jog Dial-Knob to set the value of Duration. The Duration can be set up to 99 hours 59 minutes.
- Push the MAIN button after finishing the setting of the value of Duration to store input values of the STEP and to continue the setting of the next STEP.
- Push the Jog Dial-Knob for over 1 second after completing the whole desired setting of each STEP to finalize the setting of each STEP of the selected Program no.
- If the Jog Dial-Knob is pushed shortly for once in the Set Program Mode being sustained then the set Program no. will be initiated.
- The programmed operation will not be started unless all values of the Set temperature, Rising Time, and Duration are set.
- During the programmed operation, the number of the STEP currently in operation will be displayed together with the values of current temperature and remaining time of the set Duration.
- In the Program Mode, it would be impossible to get into another mode; and to stop the operation, push the Jog Dial-Knob once shortly.
- The stored Program can be accessed easily and will enable users to start the operation conveniently by identifying the Program No. after entering into the Program Mode.

Step 0 - Temp SV : 500°C  
 Time RV : 00:05 (① Section)  
 Time MV : 01:00 (② Section)

Step 1 - Temp SV : 800°C  
 Time RV : 00:10 (③ Section)  
 Time MV : 01:00 (④ Section)

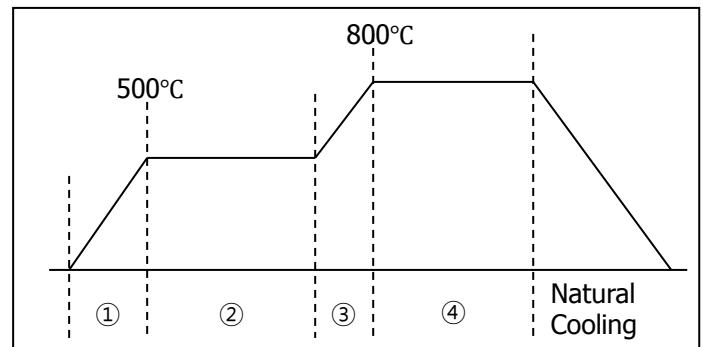- This product is different from time to reach the target temperature according to the mounting Heater capacity (Refer to specification of "Required time to be reached 800°C). If your set times in 1 and 3 were not enough to reach the target, RV value will don't start. And heating will work continuously until  $\pm 10\%$  of target temperature. When after reach the  $\pm 10\%$  of target temperature, 2 and 4 which are Time MV will start operation.

## 6.6 Key lock

- To avoid accidental changing of the set values, push the MAIN button in the menu item "set turn-off delay time".
- In this menu item can you see the temperature in the SV-line and the remaining time of delay function in the PV-line. The function of the Jog-Dial knob is deactivated.
- Press the MAIN button once more to leave the key lock. Than you are in the menu item "set speed".

## 6.7 Storage function

The **Programmable Muffle Furnace** includes a storage function for temperature and delay-timer values. That is practical if you want to work with the same parameters often. Then you need not to set them in every usage.

- For storing values you only need to hold the Jog-Dial knob pushed for two seconds. Successful storing is displayed by blinking of the SV-value for three times.
- If you saved a temperature value, "HEATING" is blinking after the device's restart.
- If you saved the delay time it is not showed at the LCD.
- Every stored parameter is erasable. Turn the value with the Jug-Dial till "0" and then hold the knob pressed for two seconds.
- If you want to store a new value it is not necessary to erase the old one.

## 7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device and pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

## 8 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> <li>✓ Check the supplied voltage is present.</li> <li>✓ Check the power connection and the power cord for proper connection and integrity.</li> <li>✓ Check if the MAIN Power Switch is "ON".</li> <li>✓ Fuse is blown.</li> </ul>
Unit Not Heating	<ul style="list-style-type: none"> <li>✓ Check if the MAIN Power Switch is "ON".</li> <li>✓ Check if the set temperature (SV) confirmed with the Jog-Dial knob (by turn the Jog-Dial knob cannot change the set temperature – see point 6.4)</li> <li>✓ Check the Safety switch is set higher (min. 40°C-50°C) than the set temperature (SV).</li> </ul>
Heating, but unit not reaching temperature or temperature is fluctuated	<ul style="list-style-type: none"> <li>✓ Check the Safety switch is set higher (min. 40°C-50°C) than the set temperature (SV).</li> <li>✓ Check the Ventilation openings are closed.</li> <li>✓ Problem in the sensor or main controller.</li> </ul>
ERROR 1	<ul style="list-style-type: none"> <li>✓ Temperature Sensor Error.</li> </ul>
ERROR 3	<ul style="list-style-type: none"> <li>✓ Door open alarm and appears when you open the door.</li> <li>✓ If Error 3 develops during door closed, check the Limit Switch is pressed properly.</li> </ul>

**If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.**

**Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.**

## 9 Technical data

	FP-03	FP-05	FP-12	FP-14	FP-27	FP-63
	FHP-03	FHP-05	FHP-12	FHP-14	FHP-27	FHP-63
<b>Chamber(cm)</b>	13×25×9	15×30×10	20×30×20	25×35×16	30×30×30	40×40×40
<b>External (cm)</b>	41x48x53	44x53x55	50x55x66	55x59x62	60x59x76	100x73x92
<b>Capacity</b>	3 Lit.	4.5 Lit.	12 Lit.	14 Lit.	27 Lit.	63 Lit.
<b>Temp. Range</b>	300 °C – 1000 °C (FHP model up to 1200°C)					
<b>Time required at 800 °C</b>	16 min		20 min		23 min	
<b>Internal</b>	Ceramic Fiber with Heating wire					
<b>External</b>	Powder Coated Steel					
<b>Controller</b>	Digital PID Controller with Jog-Dial Switch (Turn + Push)					
<b>Display</b>	Digital LCD Display with Back-Light Function					
<b>Timer</b>	99hr 59min (Continuous Function Included)					
<b>Thermocouple</b>	K-type					
<b>Heater</b>	Ceramic Fiber with Heating Wire, 4-Side Heating					
<b>Insulation</b>	Ceramic Fiber					
<b>Safety Device</b>	Over Temp. and Over Current Protector, Sensor Error Detector					
<b>Others</b>	Locking Mode (Jog-Dial Input disabled) Alarm ( Error Status and Timer-end )					
<b>Power</b>	1.4 kW	1.8 kW	3.6 kW	4.0 kW	4.6 kW	9.0 kW
<b>Packing Size (wxdxh) cm</b>	63x54x77	68x57x78	70x63x89	74x68x85	74x73x99	116x90x110
<b>Shipping Weight</b>	47 kg	54 kg	59 kg	73 kg	83 kg	225 kg
<b>Power Supply</b>	AC 230V, 1N~, 50/60Hz			AC 400 V, 3~, 50/60 Hz		

## 10 Certificate

# Certificate

for

**The quality and all features were checked by the manufacturer before the shipment.**

We grant from date of purchase

**two years guaranty.**

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

<b>Article</b>	<b>Programmable Muffle Furnace</b>
<b>Typ</b>	
<b>Serialno.</b>	
<b>Date</b>	

**witeg Labortechnik GmbH**  
Am Bildacker 16  
D-97877 Wertheim  
TEL: +49-9342-9301-0  
FAX: +49-9342-9301-77  
Email: info@witeg.de  
[www.witeg.de](http://www.witeg.de)