HB

Blockthermostat
Bedienungsanleitung
Heating Block
Operating Manual

Gilt für /written for
HB- R48, 96D

witeg Labortechnik GmbH

(English manual at page 9)

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Contents

| | |
|--|-----------|
| 1 Introduction | 10 |
| 2 Safety Instructions | 10 |
| 2.1 Description of symbols..... | 10 |
| 2.2 Instructions | 10 |
| 3 Product description | 11 |
| 3.1 Key features..... | 11 |
| 3.2 General survey | 12 |
| 3.3 Term description | 12 |
| 4 Packing contents..... | 13 |
| 4.1 Standard scope of delivery..... | 13 |
| 4.2 Additional accessories | 13 |
| 5 Positioning the device and commissioning..... | 13 |
| 6 Job description | 14 |
| 6.1 Menu structure | 14 |
| 6.3 Set temperature..... | 14 |
| 6.4 Justify | 14 |
| 6.5 Key lock | 14 |
| 6.6 Storage function..... | 15 |
| 7 Maintenance and cleaning..... | 15 |
| 8 Technical data..... | 15 |
| 9 Certificate..... | 17 |

1 Introduction

Thank you for purchasing our **HB Heating Block**. You did a good choice buying a **Wisdom** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a safe place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instructions to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid accidents caused by high temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Reparations and amendments should only be made by **witeg** authorized people. Contraventions exclude the warranty and may cause personal injury or material damages.

- Do not use this device in areas with potentially explosive atmosphere.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.

- Do not put heat-sensitive items on the device.
- Do not clean the scorching device.

- Make sure that the device can not slip away.
- To avoid heat congestions or fires make sure there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

3 Product description

3.1 Key features

The **Wisdom HB Heating Block** allows an accurate heating and a convenient and fast operation. Numerous accessories and a simple justification make this heating block to an acclaimed laboratory device in chemical, biological, medical and pharmaceutical purviews.

Key features are as follows:

- Long life by simply and solid construction.
- Simplest operating by MODE button and patented Jog-Shuttle knob.
- Turn-off delay from one minute up to over 99 hours in one minute steps adjustable.
- Signal sound after end of turn-off delay.
- Storage function for timer and temperature values.
- Illuminated LCD.
- Simple justification.
- Acrylic glass lid.
- Key lock against inadvertent changing of set speed and timer values.

3.2 General survey

3.3 Term description

SV: set value (upper line in LCD)

PV: present value, measured by the device (lower line in LCD)

Timer lamp

- Is blinking in menu items for timer modulation.
- Shines continuous in the key lock menu item when turn-off delay time runs.

Heating lamp

- Shines if the heating is activated.

Jog-Shuttle knob

- Change SV-values by turning.
- Push the knob for start and stop or probate and save set values.
- Also used for keyboard shortcut.

MODE button

- For changing between the different menu items.

Key combination

- Push the MODE button first (hold it) and in addition the Jog-Shuttle knob to reach the sub menu items.

4 Packing contents

4.1 Standard scope of delivery

| Quantity | Denomination |
|----------|--|
| 1 | HB Heating Block |
| 1 | Power supply cord |
| 1 | Acrylic glass cover |
| - | Anodized aluminium block, depending on article no. |
| 1 | Operating manual |

4.2 Additional accessories

| Order number | Denomination |
|--------------|---|
| DH.WHB510596 | BLC596 , 0,2 ml x 96 |
| DH.WHB510570 | BLC570 , 0,5 ml x 63 |
| DH.WHB511548 | BLC548 , 1,5 ml x 48 |
| DH.WHB521248 | BLT248 , Ø 12 mm x 48 |
| DH.WHB521335 | BLT335 , Ø 13 mm x 30 |
| DH.WHB521535 | BLT535 , Ø 15 mm x 30 |
| DH.WHB521628 | BLT628 , Ø 16,5 mm x 24 |
| DH.WHB521828 | BLT828 , Ø 18 mm x 24 |
| DH.WHB531515 | BLC515 , 15 ml conical x 15 |
| DH.WHB535008 | BLC008 , 50 ml conical x 8 |
| DH.WHB500101 | BLE101 , Block-extraction tong (included in every optional block) |


5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Remove perhaps soiling and package material.
5. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
6. Make sure that you follow all safety instructions.
7. Connect the device with the power outlet.
8. Switch the device on and start the process (look at job description).

6 Job description

6.1 Menu structure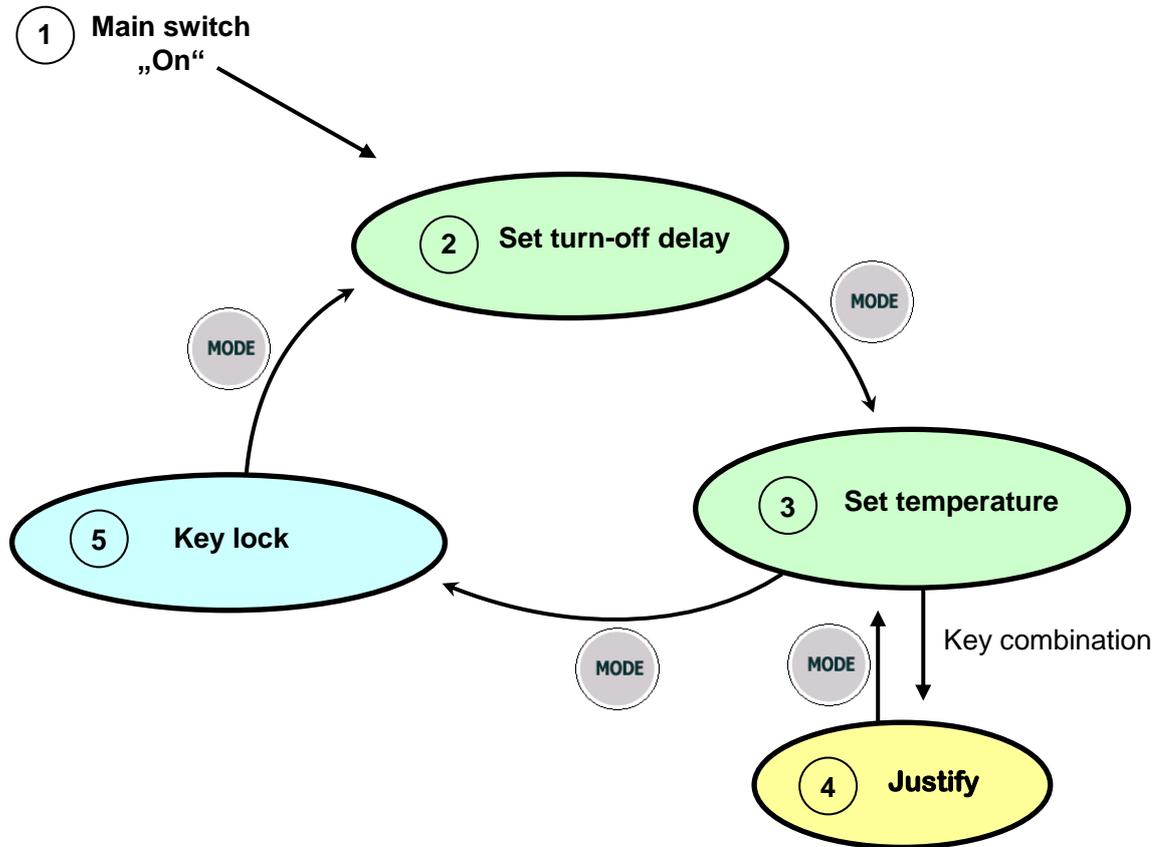

6.2 Set turn-off delay

The turn-off delay timer serves as automatically stop function. The heating will be stopped if time is up.

- Please use menu structure chart to open this menu item.
- Turn Jog-Shuttle to set timer value.
- Start and stop timer by pushing Jog-Shuttle knob.
- Value is only adjustable when timer is stopped.
- Timer lamp is blinking in this menu item.

6.3 Set temperature

- Please use menu structure chart to open this menu item.
- Turn Jog-Shuttle to set temperature value.
- Start and stop heating by pushing Jog-Shuttle knob.
- Temperature is only adjustable when heating is stopped.
- Heating lamp is blinking in this menu item.

6.4 Justify

Use the offset to compensate deviations between real and displayed temperature.

- The Offset is calculated as follows:

Temperature measured by you – showed temperature (PV) = Offset

- Please use menu structure chart to open this menu item.
- Turn Jog-Shuttle to set offset value.
- Verify offset value by pushing Jog-Shuttle knob.

6.5 Key lock

The key lock function avoids accidental change of operational parameters. Jog-Shuttle knob is deactivated.

- Please use menu structure chart to open this menu item.
- Press the MODE button to leave the key lock.

6.6 Storage function

Storage Function is very useful if you often want to work with the same parameters. Timer, temperature and offset values can be stored. Stored values will be shown when you open the according menu items and only need to be confirmed by pushing the Jog-Shuttle knob.

- Just hold Jog-Shuttle pushed for 3 seconds to save the actual value.
- Set value to "0" and hold Jog-Shuttle pushed for 3 seconds to erase stored value.
- You need not to delete a stored value to save a new one.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.


8 Technical data

| | | HB-R48 | HB-96D |
|----------------------------------|---------------|---|--------------------------------------|
| Dimension (w x d x h) (mm) | Block socked | 90x120x60 | (90x120x60) x 2 |
| | Housing | 185x230x220 | 285x250x155 |
| Capacity | | 1 | 2 |
| Heating power | | 300 W | 600 W |
| Peltier cooler | | 2 Modules | - |
| Temp. | Range | -5 °C - 95 °C | Ambient temperature +5 °C -150 °C |
| | Accuracy | ±0,1 °C at 37 °C, ±0.2 °C at 90 °C, ±0,4 °C at 135 °C | |
| | Heat up speed | ~ 5 °C/min | |
| Controller | | Digital PID controller with Jog-Shuttle knob (turn + push) | |
| | | Single Digital Fuzzy Controller | Dual Digital Fuzzy Controller |
| Display | | Illuminated LCD | |
| Timer | | 99 h, 59 min for turn-off delay | |
| Sensor | | PT100 | |
| Material | Block socked | Stainless steel | |
| | Housing | Powder coated steel | |
| | Block | Anodized aluminium | |
| | Cover | Acrylic glass | |
| Safety Device | | Over temp. and over current protector, sensor control | |
| Others | | <ul style="list-style-type: none"> - Storage function (temp., timer) - Key lock (Jog-Shuttle input disabled) - Simple justification - Signal sound (sensor error and timer-end) | |
| Power Supply | | AC 230 V, 1N~, 50/60 Hz | |

9 Certificate

Certificate

for

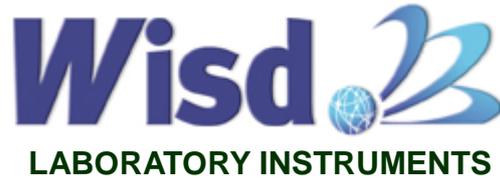The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

| | |
|------------------|-------------------------|
| Article | HB Heating Block |
| Serialno. | |
| Date | |

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de